

ADDITIONAL ADAPTER FOR PIPETTES CALIBRATION



Additional adapter for pipettes calibration is designed for gravimetric calibration and/or checking of piston pipettes. The adapter can be installed in RADWAG balances series MYA 21, XA 3Y, XA 3Y.A, XA.4Y, XA.4Y.A, AS 3Y, AS.R.

The adapter has been designed to minimize evaporation occurring in liquid handling procedures. An integral part of the adapter is an "evaporation ring" which provides high humidity rate inside the weighing chamber and the measuring vessel. Such design prevents from evaporation of liquids during weighing process.

The adapter for pipettes calibration is equipped with a specially designed weighing pan for central locating of the measuring vessel inside the weighing chamber. The whole adapter is enclosed in a glass shield equipped with a top opening for dosing liquids.

Non-centrally located opening in the top of the adapter simplifies insertion of a pipette with liquid. This solution also limits occurrence of air breeze.




The content of the adapter:

- base element;
- special weighing pan;
- measuring vessel;
- "evaporation curtain" vessel;
- glass cover with dosing opening


For balances of XA series (XA.3Y, XA.4Y, XA.4Y.A) apply pipettes calibration adapter (100 ml) only for models with max capacity higher than 100g, this is conditioned by about 65g narrower weighing range (beaker weight).

Compatible with:


- **11 ml adapter:** MYA.xx series;
- **17 ml adapter:** XA.3Y, XA.3Y.A, XA.4Y, XA.4Y.A, AS.3Y and AS.R series
- **100 ml adapter:** XA.3Y, XA.4Y (excluding XA52.3Y and XA52.4Y), XA.3Y.A, XA.4Y.A (excluding XA52.3Y.A and XA52.4Y.A), AS.3Y and AS.R series.

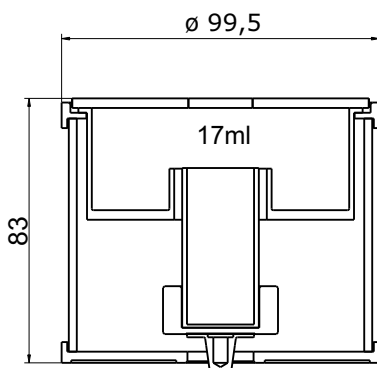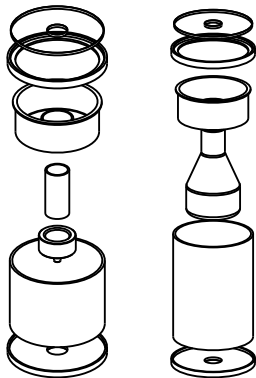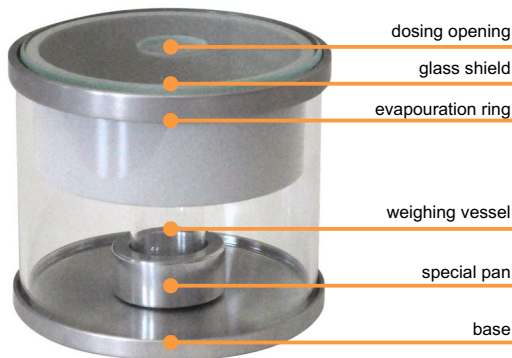

XA Series

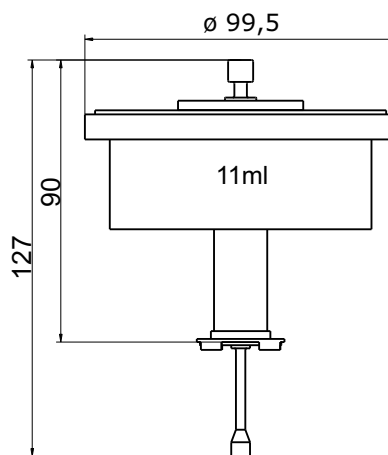

MYA Series

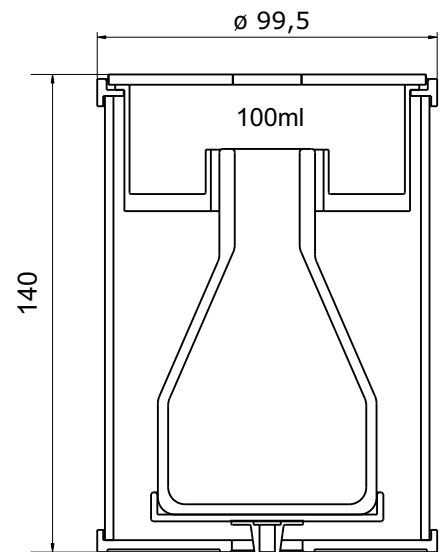

XA Series

PARTNER CORPORATION

Sos. Chitilei, Nr.16, Sector 1, Bucuresti * Tel/fax : 021-7955505 * e-mail:office@partner.com.ro * www.balante.ro